IKA® Vortex 2

Betriebsanleitung	DE	03
Ursprungssprache		
Operating instructions	EN	08
Mode d'emploi	FR	13
Руководство пользователя	RU	18
使用说明	ZH	23
取扱説明書	JA	28
사용 설명서	KO	33

Indicaciones de seguridad	ES	38
Veiligheidsinstructies	NL	39
Norme di sicurezza	IT	40
Säkerhetsanvisningar	SV	41
Sikkerhedshenvisninger	DA	42
Sikkerhetsanvisninger	NO	43
Turvallisuusohjeet	FI	44
Normas de segurança	PT	45
Wskazówki bezpieczeństwa	PL	46
Bezpečnostní pokyny	CS	47

Biztonsági utasítások	HU	48
Varnostna opozorila	SL	49
Bezpečnostné pokyny	SK	50
Ohutusjuhised	ET	51
Drošības norādes	LV	52
Nurodymai dėl saugumo	LT	53
Инструкции за безопасност	BG	54
Indicații de siguranță	RO	55
Υποδείξεις ασφάλειας	EL	56

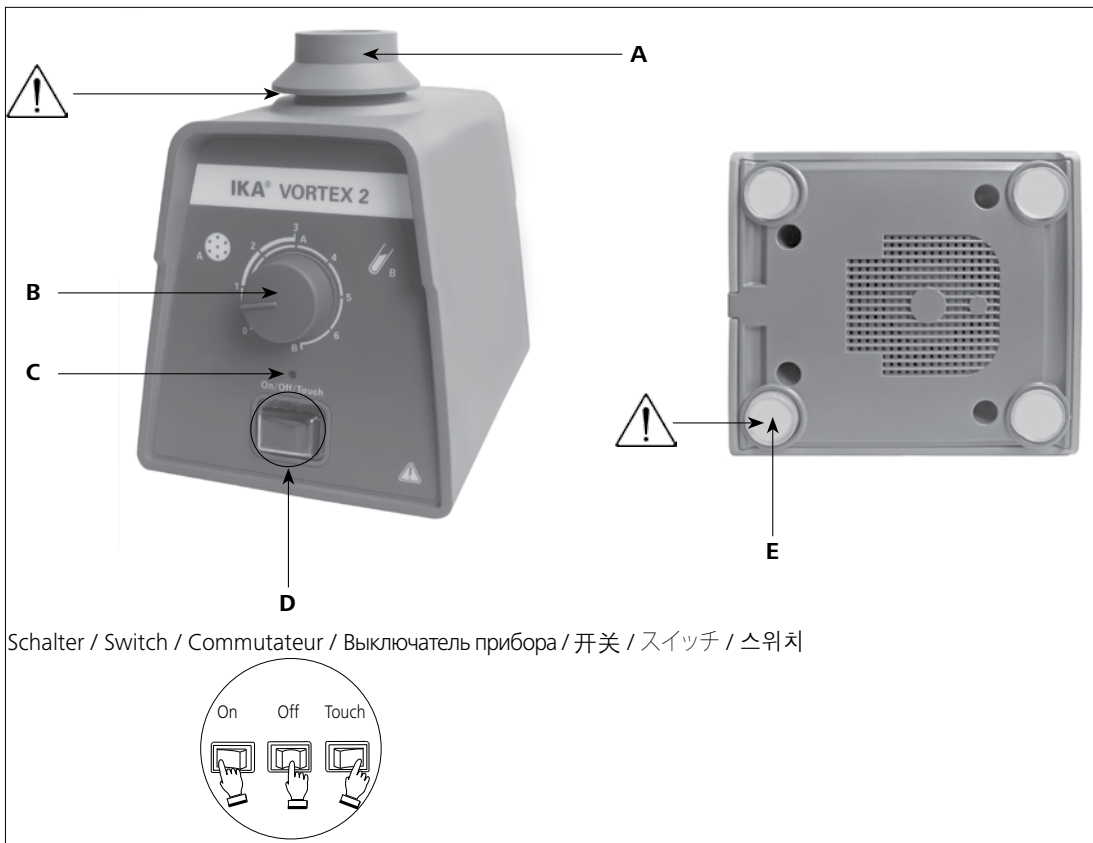
Fig. 1

Pos. Bezeichnung	Item Designation	Pos. Désignation	Пункт описание
A Standardaufsatz VG 3.1	A Standard attachment VG 3.1	A Support standard VG 3.1	A Стандартное крепление VG 3.1
B Drehknopf	B Rotating knob	B Bouton rotatif	B ручку вращения
C Kontrollleuchte	C Control light	C Lumière côntrôl	C Индикаторная лампа
D Schalter	D Switch	D Commutateur	D Выключатель прибора
E Gerätefuß	E Base	E Pied	E база устройства

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Contents

	Page
EC-Declaration of conformity	8
Explication of warning symbols	8
Safety instructions	9
Correct use	10
Unpacking	10
Commissioning	10
Accessories	11
Maintenance and Cleaning	12
Warranty	12
Technical data	12

EC-Declaration of conformity

We declare under our sole responsibility that this product is in compliance with the regulations 2006/95/EC, 2004/108/EC and 2011/65/EU and conforms to the standards or standardized documents: EN 61010-1, EN 61010-2-051 and EN 61326-1.

Explication of warning symbols

General hazard

This symbol identifies information **that is of vital importance for protecting your health and safety**. Disregarding this information may lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**.

Disregarding this information may result in damage to the instrument or to system components.

This symbol indicates information **which is important for proper use and ensuring that the operations of the instrument are performed efficiently**.

Failure to observe this information may result in inaccurate results.

Safety instructions

- **Read the operating instructions completely before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Because the options for combining products, attachments, shaking vessels, experiment and medium are nearly endless, user safety cannot be ensured simply with design requirements on the part of the product.

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise

there is a risk of:

- splashing liquids;
 - projectile parts;
 - body parts, hair, clothing and jewellery getting caught.
- Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
 - Keep the sample container as vertical as possible.
 - Only exert the minimum pressure necessary for the mixing process.
 - Use attachments for multiple samples when there are a large number of samples.
 - Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

Please pay attention to the dangerous parts of the equipment in **Fig. 1**.

The feet of the appliance must be clean and undamaged.

- If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
- Make certain that the unit is set at the lowest speed before commissioning; otherwise, the unit will begin running at the speed set in last operation. Gradually increase the speed.
- Pay attention when setting the speed to any imbalance of the attachments, shaking vessels and possible spraying of the medium to be mixed.
- Ensure that the shaker does not start to move.

Reduce the speed if:

- the medium splashes out of the vessel because the speed is too high;
- the appliance is not running smoothly;
- the appliance begins to move around because of dynamic forces;
- an error occurs.

Do not touch rotating parts during operation!

- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.

- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
- The appliance is not suitable for hand-held operation.
- Sharp edged vessels cause wear debris on the attachments.

Beware of the risk of:

- flammable materials;
 - glass breakage as a result of mechanical shaking power.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
 - **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
 - Process pathogenic materials only in closed vessels under a suitable fume hood. Please contact **IKA®** application support if you have any question.
 - Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.
 - Always disconnect the plug before fitting accessories.
 - The appliance starts up again automatically following a cut in the power supply.
 - The instrument can only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
 - The socket for the mains cord must be easily accessible.
 - The appliance may heat up when in use.
 - Never cover the ventilation slots or cooling fins on the motor or on the instrument.
 - **Warning !** This user manual needs to be consulted in all cases where it is used.
 - Socket must be earthed (protective ground contact).
 - The voltage stated on the nameplate must correspond to the mains voltage.
 - Protect the appliance and accessories from bumps and impacts.
 - Please observe the permitted speed for the attachments. Never set higher speed.
 - During operation please always keep the appliance be covered with the top attachment. Do not allow liquids to get into the appliance when changing attachments, if liquids accidentally splash into the appliance, stop using it immediately and send it back to **IKA®** for repair.
 - The machine must only be opened by trained specialists, even during repair. The instrument must be unplugged from the power supply before opening. Live parts inside the instrument may still be live for some time after unplugging from the power supply.

Covering or parts that are capable of being removed from the unit without accessory equipment have to be reattached to the unit for safe operation in order to prevent, for example, the ingress of fluids, foreign matter, etc.

Correct use

• Use

For mixing liquids:

- Touch mode for single test tubes
 - Continuous mode for single or several vessels
- Intended use: Tabletop device

• Range of use (Only for indoor use)

- Laboratories
- Pharmacies
- Schools
- Universities

This instrument is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the instrument is operated with accessories that are not supplied or recommended by the manufacturer;
- If the instrument is operated improperly or contrary to the manufacturer's specifications;
- If the instrument or the printed circuit board are modified by third parties.

Unpacking

• Unpack

- Please unpack the device carefully;
- In the case of any damage a fact report must be sent immediately (post, rail or forwarder).

• Delivery scope

- Shaker **Vortex 2**
- Standard attachment VG 3.1
- Operating instructions
- A warranty card

Commissioning


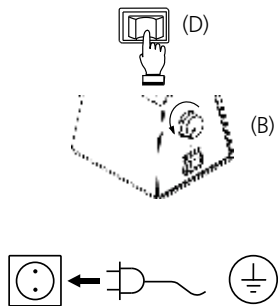
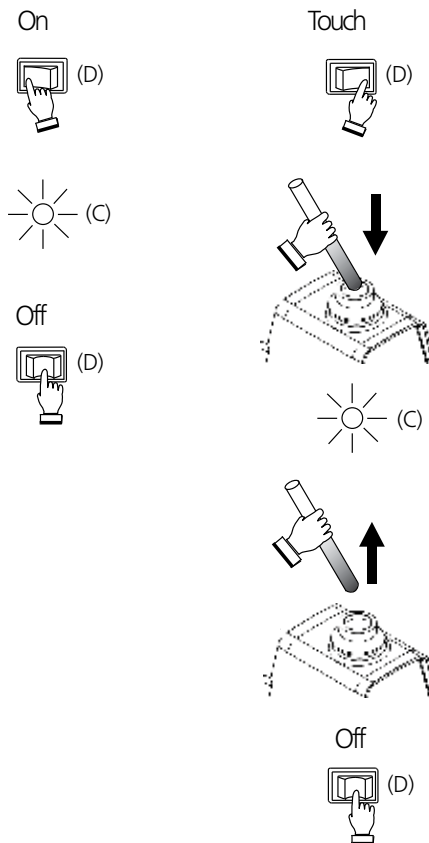
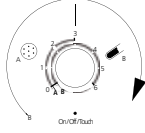
Fig. 2

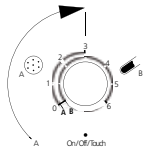
Accessories


Possible uses and permitted speed ranges of attachments:

Designation	Description	Id.-No.	Touch Mode	Continuous Mode	Range of speed Max.	Max. load (kg)
 VG 3.1 Standard attachment	For test tubes and small vessels up to diameter of 30 mm	33 412 00	×	×		0.4
 VG 3.2 One-hand attachment	88 mm-plate attachment with rubber pad	33 423 00	×	×		
 VG 3.3 Universal attachment	150 mm-plate attachment with rubber pad	33 424 00	-	×		
 VG 3.31 Test tube insert	For 54 Eppendorf tubes, for VG 3.3	33 443 00	-	×		
 VG 3.32 Test tube insert	For 18 test tubes 10 mm, for VG 3.3	33 439 00	-	×		
 VG 3.33 Test tube insert	For 12 test tubes 12 mm, for VG 3.3	33 440 00	-	×		
 VG 3.34 Test tube insert	For 8 test tubes 16 mm, for VG 3.3	33 441 00	-	×		
 VG 3.35 Test tube insert	For 8 test tubes 20 mm, for VG 3.3	33 442 00	-	×		
 VG 3.36 Erlenmeyer flask attachment	For 1 Erlenmeyer-/ round-bottomed flask 100-250 ml, for VG 3.3	33 421 00	-	×		
 VG 3.37 Microtiter plate insert	For 1 microtiter plate, for VG 3.3	33 444 00	-	×		

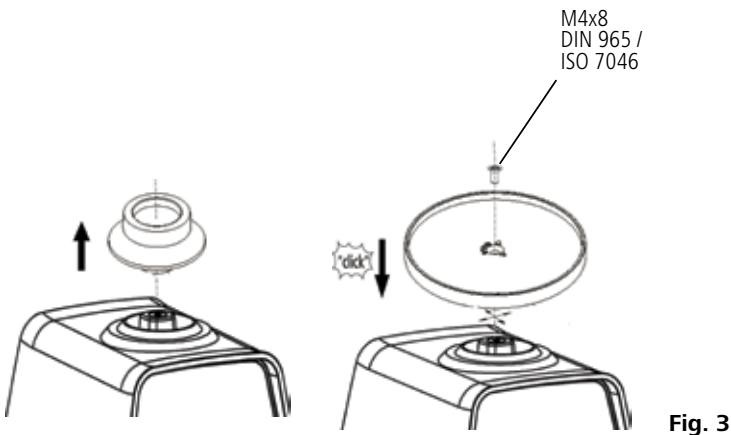Changing attachments


Using inserts

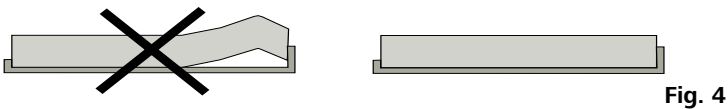Maintenance and Cleaning

The instrument is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

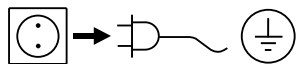Cleaning

For cleaning disconnect the main plug!

Use only cleaning agents which have been approved by **IKA®** to clean **IKA®** instruments.

Dirt	Cleaning agent
Dyes	isopropyl alcohol
Construction materials	water containing tenside/isopropyl alcohol
Cosmetics	water containing tenside/isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

For materials which are not listed, please request information from **IKA®** application support.

Wear protective gloves during cleaning the instruments.

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.

Do not allow moisture to get into the instrument when cleaning.

Before using another than the recommended method for cleaning

or decontamination, the user must ascertain with **IKA®** that this method does not destroy the instrument.

The rubber of the attachments' pads (VG 3.2 and VG 3.3) is a natural product, so talcum can discharge. Clean the rubber pad with glycerin.

Spare parts order

When ordering spare parts, please give:

- Machine type;
- Manufacturing number, see nameplate;
- Item number and designation of the spare part, see www.ika.com.

Repair

Please send instrument in for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For this you should request the "Decontamination Clearance Certificate" from **IKA®**, or use the download printout of it from the **IKA®** website www.ika.com.

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Warranty

In accordance with **IKA®** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

Technical data

Operating voltage / Design frequency	V/Hz	220...240 ±10% / 50/60
or	V/Hz	115±10% / 60
or	V/Hz	100±10% / 50/60
Input power	W	60
Speed	rpm	(750 - 2750) ± 250, at 50 Hz, infinitely adjustable (750 - 3300) ± 250, at 60 Hz, infinitely adjustable analog (infinitely) rotating knob in front of the device
Speed adjustment		rotating knob in front of the device
Speed display		Scale 0 - 6
Drive		Shaded pole motor
Agitation stroke	mm	4
Shaking motion		horizontal, approximately circular
Perm. duration of operation	%	100
Perm. ambient temperature	°C	+5 to +40
Perm. relative humidity	%	80
Protection type (acc. to DIN EN 60529)		IP 21 (IP 21 can only be achieved when fitted with top attachment)
Protection class		I (protective earth)
Overvoltage category		II
Contamination level		2
Operation at a terrestrial altitude	m	max. 2000
Dimensions total (W x D x H) (without attachment)	mm	120 x 138 x 140
Weight without supported load	kg	3.9
Max. supported load (incl. attachment)	kg	0.4

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IKA® - Werke GmbH & Co.KG